


----- clip, laminate, keep with freezer -----

### Initial Setup

1. Unpack freezer unit and included components, ensure all packaging material has been removed from around the fan inlet and outlet holes.
2. Remove orange Transport Stop and replace with the white Protective Cap before powering on the freezer.
3. Plug AC power cord into the main power socket at the back of the unit and then into an electrical outlet
4. Turn the AC Power Switch to the "On" position.
5. Notice: Ensure the Access Port stopper or sealant is in place. Remove factory tape at the port stopper prior to initial use.
6. Turn on battery power. The Battery Power Switch is located within the Recessed Electrical Panel on the rear of the freezer.
7. Ensure setpoint is within desired range (default is -80°C) or change setpoint to user-preferred temperature according to the control functions outlined below.

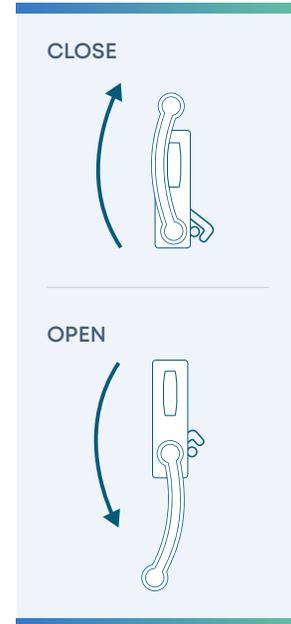

**NOTICE:**

Before engaging door latch, ensure handle is fully rotated down to the stop point in order to reset latching mechanism. To engage door latch when closing, push out door IN before rotating door handle to UP position.

### Guide to Basic GUI Functions

<b>ON / OFF</b>	Switch the main power on, located at the back of the unit above the AC cord.
<b>Battery Power</b>	Switch on the battery power, located within the Recessed Electrical Panel on the rear of the freezer. This is essential if monitoring is required in the event of a power failure.
<b>Change Setpoint</b>	Tap Home > Setup > Setup > Enter PIN > Enter > Change > Setpoint > +/- > Save
<b>Silence Alarm</b>	Tap Home > Setup > Silence
<b>Event Log</b>	Tap Home > Setup > Event Log > Prev. Page, Next Page
<b>Chart History</b>	Tap Home > Setup > Chart History

### Guide to Advanced GUI Functions

<b>Alarm Set-up</b>	Tap Home > Setup > Setup > Enter PIN > Enter > Alarms > xxx Delay > +/- > Save
<b>Date and Time</b>	Tap Home > Setup > PIN > Date & Time > +/- > Save
<b>Change PIN</b>	Tap Home > Setup > Setup > Enter PIN > Enter > PIN > Enter New PIN > Enter > Re-enter New PIN > Enter
<b>Calibrate RTDs</b>	Tap Home > Setup > About > Service > RTD Calibr. > RTD Select > +/- > Calib.

6000 Poston Road, Athens, Ohio 45701, USA  
t: 740.274.7900 / 855.274.7900 f: 740.274.7901 www.stirlingultracold.com

©2022 Stirling Ultracold, a part of BioLife Solutions.

Global Cooling technology is manufactured under U.S. and International patents. Stirling Ultracold is a trademark of Global Cooling, Inc. Specifications subject to change without notice. Refer to [stirlingultracold.com](http://stirlingultracold.com) for the latest specifications.

WE  
KNOW  
COLD™