


SPECIFICATIONS

→ Application, rating and electric data

Application	Storage of general (non-flammable) laboratory materials
Storage Volume	25 liters (0.9 cu.ft.)
Storage Capacity	18 standard 2" boxes
Temperature Range	-86°C to -20°C @ 32°C (90°F) ambient, uniformity +/- 3°C at -80°C top to bottom, adjustable in 1°C increments. Default temperature setting is -80°C. Preset options are -86°C, -40°C and -20°C.
Electric Power	100V to 240VAC at either 50 or 60Hz or 12V DC from mobile source
Maximum Power (Current)	280 watts (4 amps @120VAC, 2 amps @240VAC, 15 amps @ 12V DC)
Auto-Voltage Capability	100V to 240V, 50 or 60 Hz
Electric Supply Rating	15 amp or greater grounded circuit
Power Cord Options Available	Multiple options are available. Consult with your sales representative
	
Certification/Agency Listing	CE, cTUVus
Noise	Advanced noise abatement, <45 dB(A) at 1 meter
Indoor/Outdoor Use	Indoor use only
Environmental Conditions	Non-corrosive, non-flammable, non-explosive, indoor use, altitude up to 2,000m, temperature 5°C to 40°C (41°F to 104°F), maximum relative humidity 80% f or temperature up to 31°C decreasing linearly to 50% at 40°C

→ Refrigeration system

Cooling Engine	Helium charged free-piston Stirling engine with continuous modulation
Heat Transport System	Gravity driven thermosiphon
Refrigerant	R-170 (Ethane) 10–12 grams
Evaporator	Cold wall (inner liner)
Heat Rejection	Finned heat exchanger with forced air cooling Air inlet: Right front side of unit Air outlet: Below control panel of front mechanical compartment
Defrost Method	Manual

→ Controller


Controller Type	Microprocessor controls
Security	Lockable lid
Warm and Cold Alarms	Setpoint ± 10°C
Control Sensor	One RTD (PT100 Class A)
Dry Contacts	Optional
Internet Connectivity	Optional SenseAnywhere wireless temperature monitoring and logging

→ **Performance data**

Steady State Energy Use at 25°C Ambient	2.8 kWh/day (average power 118 watts) at -80°C (Empty Cabinet)
Pull-Down from 25°C Ambient	4 hours to -80°C (Empty Cabinet)
Recovery from 1 minute lid opening	20 minutes to -80°C (Empty Cabinet)
Warm-up Profile	30 minutes to -60°C (Empty Cabinet) 70 minutes to -40°C (Empty Cabinet)
Heat Dissipation	403 BTU/h (load to HVAC) (Empty Cabinet)

→ **Dimensions and construction**

Interior (L x W x D)	332 x 221 x 340 mm / (13.1 x 8.7 x 13.4 in.)
Exterior (L x W x D)	692.5 x 350 x 460 mm / (27.3 x 13.8 x 18.1 in.)
Net Weight, Empty	21 kg (46 lbs.), nominal
Shipping (L x W x D)	762 x 457 x 559 mm / (30 x 18 x 22 in.)
Shipping Weight	24 kg (53 lbs.)
Insulation	High performance vacuum insulated panels and polyurethane foam
Gasket Heater	Continuous operation when the power cord is connected
Options	Plastic Utility Bin, AC Power Cords, Mobile Power DC Cord, Sub-Lid, Certificate of Calibration, Temperature Mapping Validation, Remote Alarm Contacts

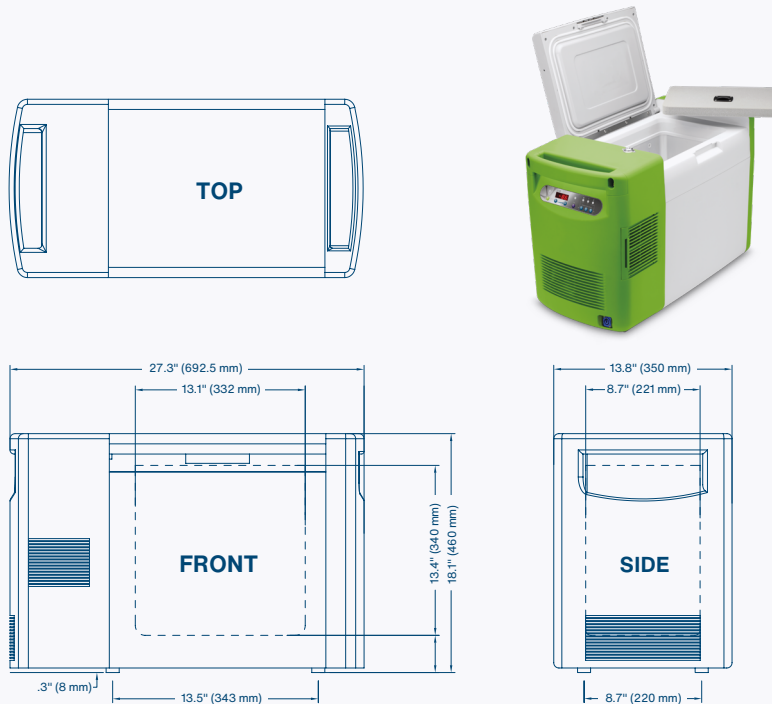

6000 Poston Road
Athens, OH 45701 USA
T +1.740.274.7900 / 855.274.7900
F +1.740.274.7901
E info@stirlingultracold.com
BioLifeSolutions.com

2023© Stirling Ultracold,
a part of BioLife Solutions.
All rights reserved.

Global Cooling technology is manufactured under U.S. and International patents. Stirling Ultracold is a trademark of Global Cooling, Inc. Specifications subject to change without notice.

Storage Solutions

SAL-SPC-00004-D